



MEDISCA®
YOUR TRUSTED PARTNER IN COMPOUNDING

RÉFÉRENCE DU PRODUIT : 6838-01



PIÈCES DE RECHANGE, ProFiller 3600 (TAILLE #000)

DESCRIPTION : Les pièces de rechange* pour une taille de capsule comprennent : plateau à capsules / jeu de tôles de corps (A, B, C) / plaque d'écartement / plateau d'orientation / casier à capsules

INFORMATIONS COMPLÉMENTAIRES :

*Un pilon peut être nécessaire. Veuillez communiquer à votre représentant commercial Medisca le numéro de série FB (situé à l'arrière de la base de remplissage) ainsi que les tailles de capsules actuellement utilisées sur votre machine à capsules afin de vous assurer que toutes les pièces nécessaires au remplissage correct des capsules sont fournies.

Produits associés :

6837-01 : Poussoir, 60 broches, taille #000

Installation des pièces de rechange :



**No Tools
Required**

ProFiller 3600 system's patented design requires no tools or adjustments to change capsule size.



1. Lift out Orienter Tray.




2. Slide new Orienter Tray in place.




3. Unscrew 4 thumbscrews from corners of Filler Base.



4. Remove Body Sheet Set.

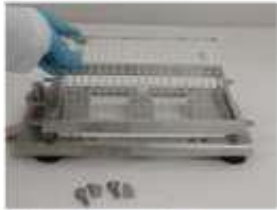

5. Switch Spacer Plate.

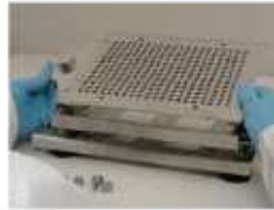

6. Position new size Body Sheet Set on Filler Base.



7. Secure with four thumb screws.